



Procedure for using the Cozap™ Coomassie® Blue Destaining Pads (0070-3717)

1. Place one of the Cozap pads in a tray with your destaining solution (25% Methanol) and gel(s).
2. Cozap pads will absorb Coomassie Blue, leaving only the protein features.
3. Once the gel is destained, remove it from the tray.

Notes:

- Cozap pads can be used for several gels, until it turns a deep blue color.
- When replacing old Cozap pads with new, it is not necessary to change the destaining solution.
- Keep Cozap pads stored at room temperature.

Cozap Coomassie Blue Destaining Pads

0070-3717 5 per pack

Collodial Coomassie Blue Stain

0080-0216 1 L

0080-0017 100 ml

Coomassie® is a registered trademark of Imperial Chemical Industries, PLC
Genomic Solutions™ and Investigator™ are trademarks of Genomic Solutions,
Inc.
Cozap™ is a trademark of Sialomed, Inc.

©2002 Genomic Solutions Inc.

